

# Protein Detection Kit for Electrophoresis Gels

---

**ENZ0003**

***Peace of mind (2 in 1)<sup>TM</sup>***

## Instructions for use

---

- At the end of the electrophoresis run, remove your gel in the usual way;
- Remove the lid from the ***Peace of mind (2 in 1)<sup>TM</sup>*** tray and place the gel (carefully) over the black layer in the bottom of the tray;
- Pour entire content of one bottle of **Stain/Destain Solution** into the tray with the gel;
- Ensure the gel is flat and completely covered with solution;
- Replace the lid (tightly) on the tray and place the sealed ***Peace of mind (2 in 1)<sup>TM</sup>*** tray on a shaking platform;
- Shake gently for **12 - 15 hrs.**

### *Step 1*

- Open the tray and remove the gel from destaining solution/tray;
- Store the gel in deionised water or dry in gel-dryer for permanent storage.

### *Step 2*

---

**NB:** Please follow the instructions carefully and precisely to obtain best results. Correct timing and use of designated solutions are essential for good results.

#### **WARNING**

**All solutions contain acetic acid and methanol and should be handled with care and disposed of in an appropriate manner.**

---

For more information please contact: Enzolve Technologies Limited  
Belfield Innovation Park, UCD  
Belfield, Dublin 4, Ireland  
Tel. (+353 1) 716 3633  
Fax (+353 1) 716 3632  
E-mail: [info@enzolve.com](mailto:info@enzolve.com)

***Enzolve***

---