

Protein Detection Kit for Electrophoresis Gels

ENZ0003

Peace of mind (2 in 1)TM

Instructions for use

- At the end of the electrophoresis run, remove your gel in the usual way;
- Remove the lid from the ***Peace of mind (2 in 1)TM*** tray and place the gel (carefully) over the black layer in the bottom of the tray;
- Pour entire content of one bottle of **Stain/Destain Solution** into the tray with the gel;
- Ensure the gel is flat and completely covered with solution;
- Replace the lid (tightly) on the tray and place the sealed ***Peace of mind (2 in 1)TM*** tray on a shaking platform;
- Shake gently for **12 - 15 hrs.**

Step 1

- Open the tray and remove the gel from destaining solution/tray;
- Store the gel in deionised water or dry in gel-dryer for permanent storage.

Step 2

NB: Please follow the instructions carefully and precisely to obtain best results. Correct timing and use of designated solutions are essential for good results.

WARNING

All solutions contain acetic acid and methanol and should be handled with care and disposed of in an appropriate manner.

For more information please contact: Enzolve Technologies Limited
Belfield Innovation Park, UCD
Belfield, Dublin 4, Ireland
Tel. (+353 1) 716 3633
Fax (+353 1) 716 3632
E-mail: info@enzolve.com

Enzolve
